

Determination of Enzyme Kinetic Parameters in Human Liver Microsomes or Supersomes™

1. Objective:

The determination of enzyme kinetic parameters in human liver microsomes or Supersomes™ is designed to provide information about the involvement of the different human CYP isoenzymes in the metabolic fate of a test compound.

2. Introduction

Liver microsomes are subcellular fractions (mainly endoplasmatic reticulum) containing many drug-metabolizing enzymes, e.g. cytochrome P450s (CYPs), flavin-monoxygenases, carboxylesterases, and epoxide hydrolase. Therefore they are widely used as an *in vitro* model system in order to investigate the metabolic fate of xenobiotics. The most prominent group of drug metabolizing enzymes is the super family of cytochrome P450s (CYPs) These haem-containing enzymes play a key role in the metabolism (mainly oxidation) of a variety of chemically diverse compounds including food compounds, pharmaceutical agents, carcinogens, and environmental pollutants.

Human liver microsomes contain the following CYP isoenzymes involved in drug metabolism: CYP1A2, 2A6, 2B6, 2C8, 2C9, 2C19, 2D6, 2E1, and 3A4. Of these isoenzymes CYP3A4 plays a major role in metabolism of xenobiotics as it is the most abundant CYP in human liver

(approx. 28 %) and it is involved in metabolism of more than 50 % of all pharmaceuticals applied in present-day medication.

Supersomes™ (human recombinant enzymes) are microsomes prepared from insect cells infected by baculovirus and containing cDNA of a single human CYP

For more information please contact us!

isoenzyme. Therefore Supersomes™ have the advantage of unique specification by expressing enzyme activity of one single CYP isoform.

The two most important enzyme kinetic parameters are:

- K_m Michaelis-Menten constant (concentration at which 50 % of the maximal velocity is observed)
- V_{max} maximum reaction velocity of enzyme

3. Short summary: Enzyme Kinetic Parameters Procedure

a) Human Liver Microsomes

- Preparation of buffers and stock solutions of test compound (usually in DMSO or water)
- Incubation of reaction mix including human liver microsomes, test compound (multiple concentrations, usually 8 to 10), and NADPH for a single time points, e.g. 60 minutes.
- End of reactions by addition of stop reagent, sample preparation
- Determination of metabolite formation and loss of parent compound (compared to zero time point control and/or no NADPH-control) using LC-MS methods.
- Results: Determination of turnover rates and calculation of enzyme kinetic parameters (K_m , V_{max}).

For more information please contact us!

b) Supersomes™

- Preparation of buffers and stock solutions of test compound (usually in DMSO or water)
- Incubation of reaction mix including Supersomes™ (human recombinant enzymes), test compound (multiple concentrations, usually 8 to 10), and NADPH for a single time points, e.g. 60 minutes.
- End of reactions by addition of stop reagent, sample preparation
- Determination of metabolite formation and loss of parent compound (compared to zero time point control and/or no NADPH-control) using LC-MS methods.
- Results: Determination of turnover rates and calculation of enzyme kinetic parameters (K_m , V_{max}).

For more information please contact us!

4. Marker Reactions

The following marker reactions and marker substrates can be employed as quality criteria of the metabolic capability of the microsomes or the Supersomes™:

CYP	Marker Substrate	Marker Reaction
1A2	7-ethoxyresorufin	7-ethoxyresorufin O-deethylation
2A6	coumarin	coumarin 7 - hydroxylation
2B6	S-mephenytoin	S-mephenytoin N-demethylation
2C8	paclitaxel	paclitaxel 6a-hydroxylation
2C9	diclofenac	diclofenac 4'-hydroxylation
2C19	S-mephenytoin	S-mephenytoin 4'-hydroxylation
2D6	bufuralol	bufuralol 1'-hydroxylation
2E1	chlorzoxazone	chlorzoxazone 6-hydroxylation
3A4	testosterone	testosterone 6β-hydroxylation

For more information please contact us!